













Handling :

Any handling of weights and masses shall be in accordance with adapted hygiene rules.

Avoid :

- taking weights of classes E and F with bare hands
- exposing the weights to dust and dirt
- storing weights in a humid atmosphere
- submitting weights to a strong magnetic fields
- pulling or pushing weights on a surface

ACCESSORIES FOR STANDARD WEIGHTS AND MASSES OF CLASSES E AND F

	<=100 g	200 to 500 g	500 g to 2 kg	5 to 50 kg	>50 kg
Non-magnetic tweezer 12,5 cm					
Non-magnetic tweezer 20 cm					
Weight forks and handles					
Gloves					
Soft cloth					
Cleaning set					

Maintenance :

No maintenance is required.

Be carefull : Any intervention that could change the value of the mass should be avoided, for example :

- polishing
- repainting
- opening the adjusting cavity

Storage :

It is recommended to store standard weights and masses in their original boxes. Weights and masses without box shall be stored in appropriate places.



Environment	Storage of weights of classes E and F	Storage of weights of class M
Storage temperature	20°C ± 5°C	20°C ± 10°C
Humidity	50% ±25%	
Dust	Off - Except bad weather	

R-COM.IN.SC.1401 04

ZWIEBEL SAS - SAINT JEAN SAVERNE - BP 50002 - F-67701 SAVERNE CEDEX

Tel. : +33 (0)3.88.71.53.10 - E-mail : commercial@zwibel.fr - www.zwibel.fr